





Parts List							DESCRIPTION
ITEM	QTY	PART NUMBER	3 FT.	4 FT.	5 FT.	6 FT.	
1	2	A-425-7373					Lateral Biosafety Cabinet Latch
2	8	A-500-395					Basestand Leg Bushing
3	2	B-500-386					Basestand End Panel Weldment
4	4	B-500-390					Basestand Leg Extension
5	1	B-500-391-TAB	-01	-02	-03	-04	Cross Member Panel Weldment 3'
6	2	B-500-393					Front Mounting Bracket
7	2	B-500-394					Rear Mounting Bracket
8	4	X-980443					LEG LEVELER FOR 2" SQ TUBING
9	12	X-980493					Acorn Nut 1/4-20 (Low Crown) SST
10	8	X-980494-03					Acorn Nut 3/8-16 SST
11	8	X-980495					Nutsert 1/4-20
12	2	X-980659-02					Bolt 3/8-16 X 3/4 SST
13	20	X-980659-12					Bolt HHCS 1/4-20 X 5/8 SST
14	10	X-980781-07					Lock Washer 3/8 Internal Tooth SST
15	18	X-980783-07					Washer 3/8 SAE Nickel
16	32	X-980794-02					Washer 1/4 SST
17	24	X-980815-01					Lock Washer 1/4 Internal Tooth SST
18	8	X-980929-01					Bolt 3/8-16 X 2-3/4 SST

REV	ECO	DESCRIPTION	DATE	DFTM	CHKD
E	13582	UPDATED HARDWARE	11/2/2018	TH	DSH

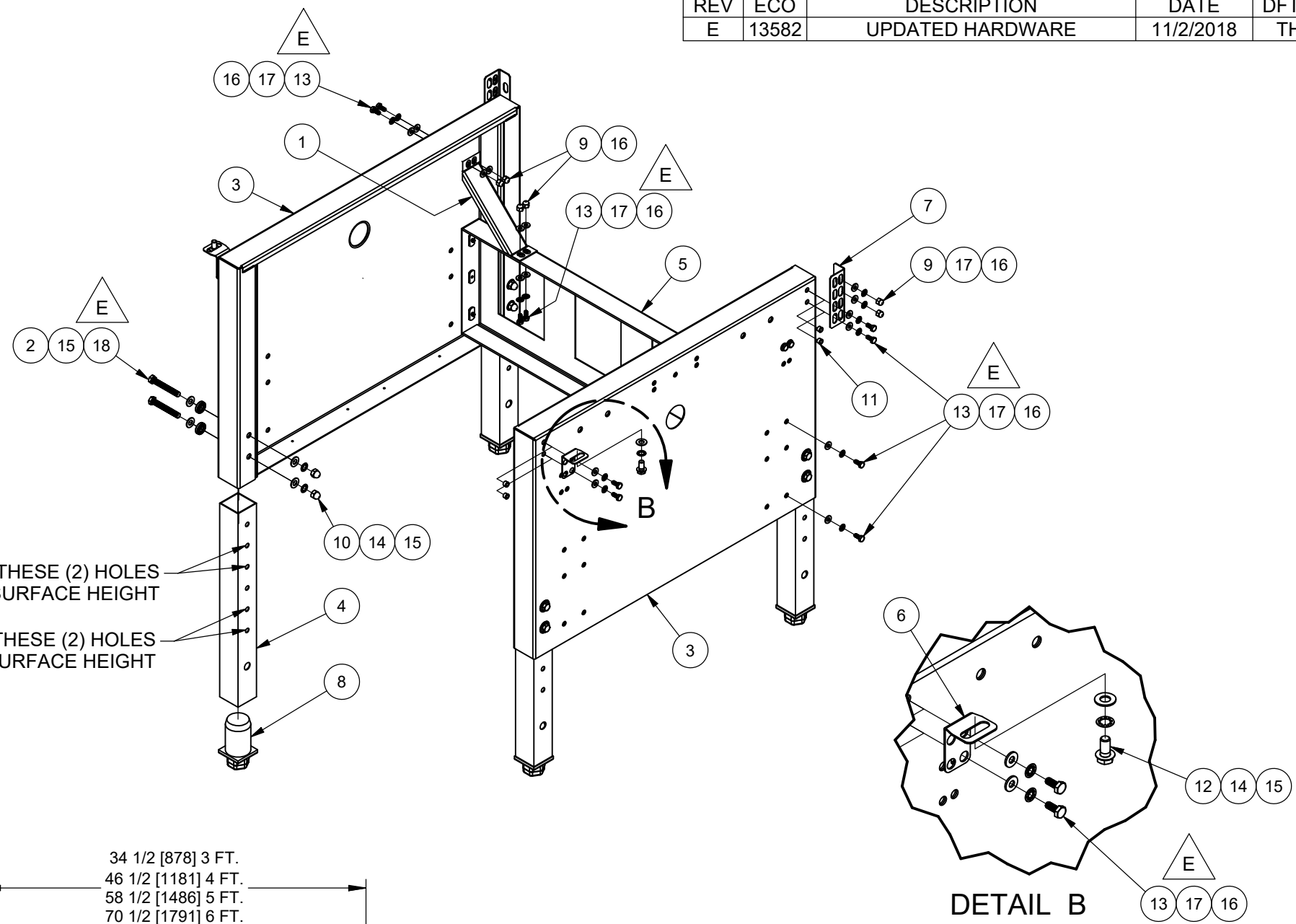


BOLT THRU THESE (2) HOLES
FOR 36" WORK SURFACE HEIGHT

BOLT THRU THESE (2) HOLES
FOR 30" WORK SURFACE HEIGHT

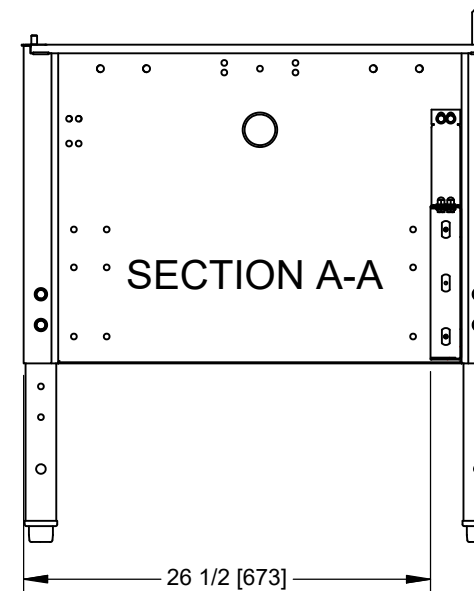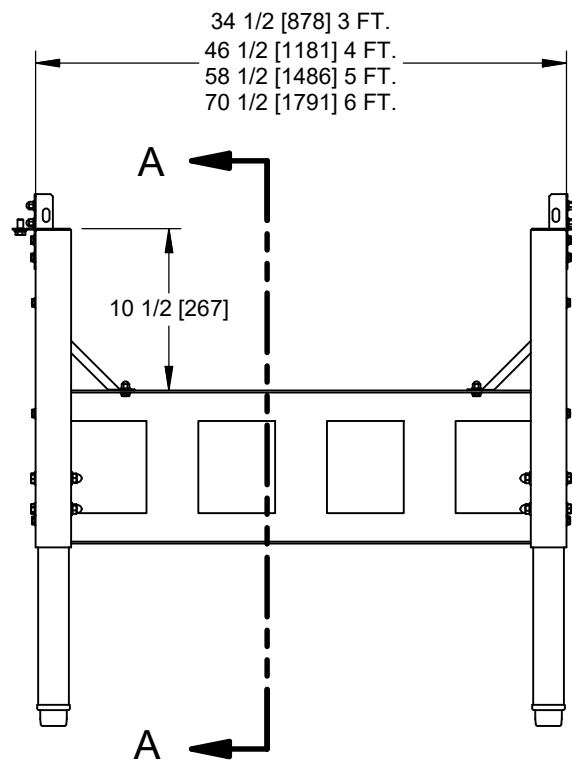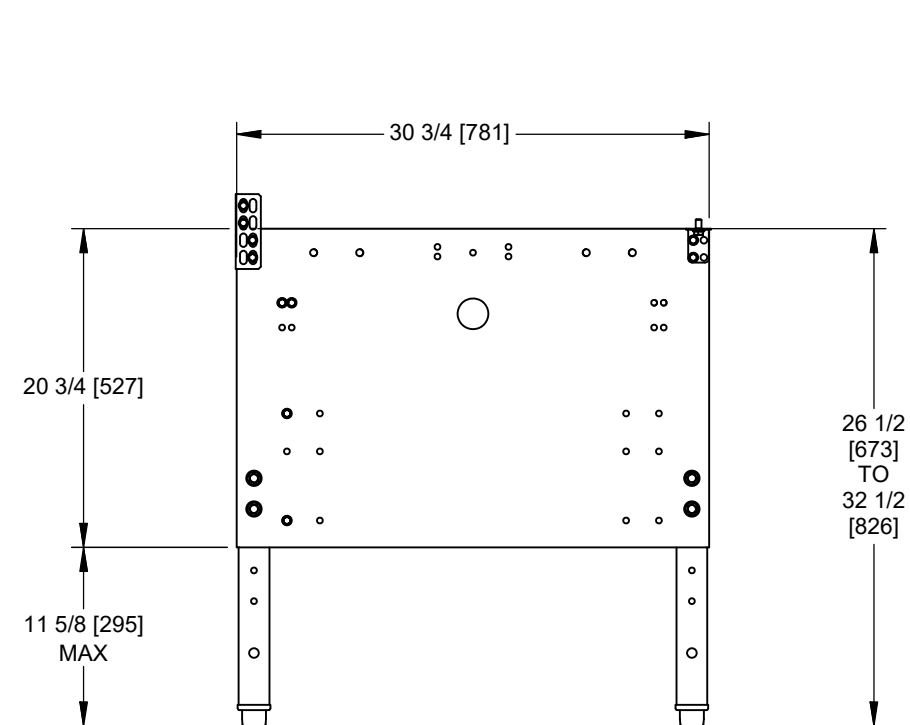

NU-475, 477, 480, 540, 543, 545, 560, 565
NU-500-100
NU-500-101
NU-500-102
NU-500-103

					
TITLE ASSEMBLY 31", TELESCOPING BASESTAND W/ LL					
DFTM LS	DATE 3/11/2014	CHKD BP	SHEET 1	OF 1	
DRAWING NUMBER BCD-16633			E		